ITCL-01

Method based on the observation of the appearance of a flocculation process, to allow measure of the process known as **test of coagulation time** of "Standard" **milk substrate**, after the addition of enzymatic products to the milk, used as coagulation agents. A known volume of enzymatic solution is mixed under fixed speed with a known quantity of milk substract, at constant and accuracy controlled temperature, and measuring the time elapsed until the observation when the flocculation process starts. The apparatus is composed by a precision thermostatic water bath, able to work at  $30^{\circ}\text{C} \pm 0,05^{\circ}\text{C}$ , with internal circulation pump, with digital temperature control; a driving system to rotate 4 test tubes at 3RPM with an inclination of  $30^{\circ}$  of the liquid level. Equipped with an electric white light of 100 Watts, to light the test tubes perfectly. To increase the contrast, there is a black surface at the rear of the test tubes. Sturdy metallic painted housing, all parts in direct contact with the water are of stainless steel.

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